



Item No.: BT3000-MR

Platform Installation:

Your Orbi-Shaker comes complete with a standard - rubber mat platform for holding flat vessels at low to medium speeds. In order to install the MAGic Clamp platform, you must follow the steps below:

- Remove the rubber mat from the shaker
- **Unscrew the 4 central screws from the attached platform: DO NOT REMOVE THE PLATFORM. (save screws for later use)**
- Remove the black mat from the MAGic Clamp platform and place the platform on the shaker so that the four center screw holes line up with the 4 holes in the shaking mechanism.
- Use the recently removed screws to screw the MAGic Clamp platform to the four center screw holes.

Flask Clamp Installation and Alignment

The MAGic Clamp Platform is capable of supporting 2 x 2000ml, 5 x 1000ml, 9x500ml, 14 x 250ml flasks, 24 x 125ml or various combinations of these flask sizes.

To attach a flask clamp, simply place on the platform in the desired location. When attaching flask clamps for the maximum capacity of the platform, please follow the diagrams below for guidance:

When attaching a mixture of different flask sizes, ALWAYS ensure a balanced load..

The MAGic Clamp platform is also capable of supporting two of the H1000-MR-1550, tube rack for 15 and 50ml tubes. When installing this tube rack simply place on the platform in the desired location .

