

Deactivate Old Sensor

1

Unplug Old Sensor

2

Plug in New Sensor

3




Add New Sensor to Account

4


To Deactivate a Sensor

1. On main screen, tap on devices.
2. Select Device to be deactivated.
3. Tap on Settings.
4. Tap Edit button.
5. Tap Deactivate, and confirm you wish to make that change.

Note: "Device" in the case of Ethernet refers to the removable sensor probe.

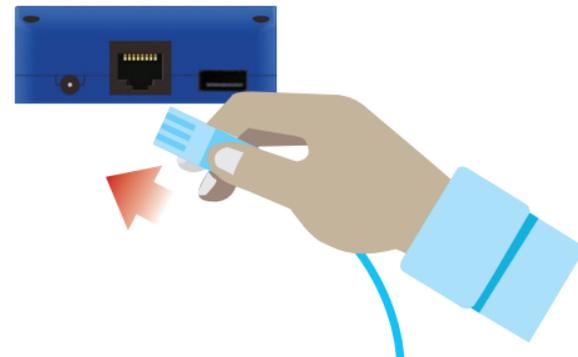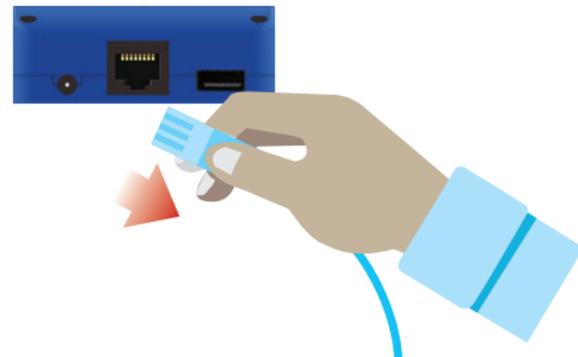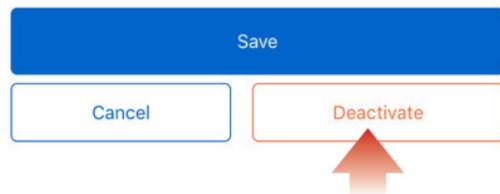

Open TraceableLIVE app on mobile or desktop and Log in to account:

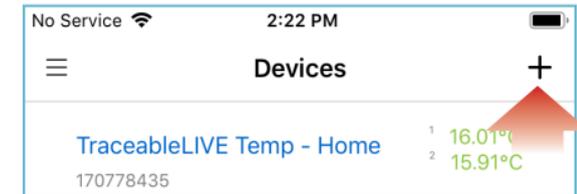

Tap the **+** icon to add a device.

- Permission must be given to TraceableLIVE app to access your mobile device's camera.
- Using your mobile device's camera hover over QR code on back of unit (see right).
- App will automatically fill in device information.



Note: QR code only works with a mobile device.

Optional: Maintain continuous data history, select a predecessor device, **see Step 5.**


Predecessor Settings

5

To maintain continuous data history from previous sensor, select a predecessor device to link data from another unit which has become inactive.

To carryover data from old sensor

1. On main screen, tap on devices.
2. Select Device to be deactivated.
3. Tap on Settings.
4. Tap Edit button.
5. Tap on drop-down menu and select device, transferring the legacy data to the new sensor.

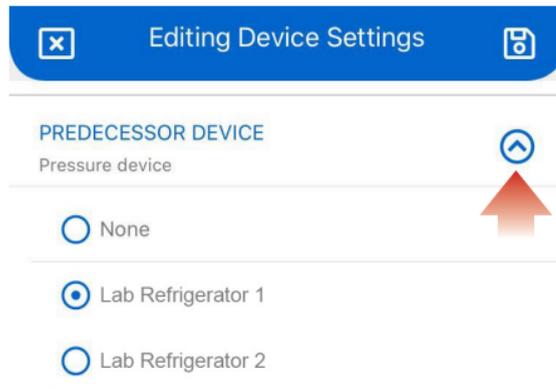

Having trouble?
Please give us a call
281-482-1714 or email
support@traceable.com

TraceableLIVE® QUICK START SENSOR GUIDE



**REPLACEMENT
SENSOR**